# Schnellverascher ULTRA X 052

## Rapid incinerator ULTRA X 052

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**This unique rapid incinerator was developed as a budget-priced alternative to kilns. It allows you to incinerate and ash single samples quickly and efficiently.**

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### ULTRA X 052 Rapid incinerator

The operation of the device is very easy. After placing the sample within the sample bowl in the preheated incineration chamber, the radiator is swung into position to start the ashing process.

The maximum temperature of approx. 610 °C can be pre-selected by in the menu. 2 heating stages are available with 2 different powers and times.

Depending on the matter of the material to be incinerated, the typical ashing time for a 5 g sample generally is between 8 and 15 minutes.

The ULTRA X 052 incinerator comes with an electronic with LCD display and 4 switches for the menu and the times and power settings.

The heating time is shown in the display and stage 2 follows automatically when stage 1 is ready. At the end of the process the heater is switched off.

A light and a signal shows when the device is in use.

### Use

This infrared rapid incinerator is used particularly in conjunction with ULTRA X laboratory devices to determine the incineration residue in sewage sludge after the determination of dry matter. The UX 052 is useful in every laboratory where a fast incineration of single samples of organic materials is needed.

To determine the dry matter and residue on ignition in sewage sludges we offer the moisture tester UX 2011 G in connection with the incinerator UX 052.

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Note:

Because noxious and/or malodorous fumes are likely to be emitted during the calcination or ashing process, it is recommended that you use the ULTRA X rapid incinerator in a protection cabinet.

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**Technical data of the ULTRA X 052**

Power supply	230 Volt, 50-60 Hz
Other voltages with separate transformer	
Radiator power:	375 Watt
Temperature range	up to approx. 610 °C
Temperature setting	electronically, 2 steps possible
Time setting	electronically, 2 steps possible up to 99 min. per stage
Sample volume	max. 60 cm <sup>3</sup> (quartzglass pan) max. 130 cm <sup>3</sup> (aluminium pan)
Max. bowl size	75 mm diameter x 30 mm height
Dimensions	Height 180 mm Width 190 mm Depth 350 mm
Weight	approx. 4.0 kg

Subject to technical changes.

Accessories available upon request:

Item	Order No.
Quartz glass bowls	0015080
Aluminium foils 0,03, Ø 130mm	1011390
Foil press to shape the aluminium foils	1230000

Our product range:

Infrared moisture testers for different usage, sample dryers (without balance), incinerators.