
Schnellverascher ULTRA X 052

Rapid incinerator ULTRA X 052

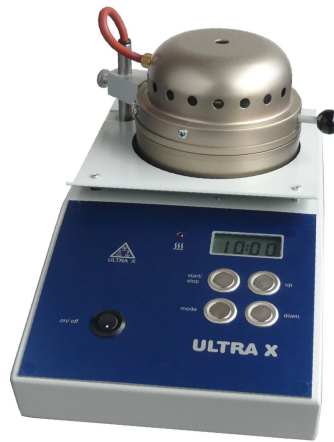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This unique rapid incinerator was developed as a cost-effective alternative to kilns. It allows you to incinerate and ash single samples quickly and efficiently.

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ULTRA X 052 Rapid incinerator

The Operation of the device is very easy. After placing the sample in the sample bowl in the preheated incineration chamber, the radiator is swung into position to start the ashing process.

The maximum temperature of approx. 610 °C can be preselected in the menu. Two heating stages are available with 2 different powers and times.

Depending on the type of material to be incinerated, the typical ashing time for a 5 g sample generally is between 8 and 15 minutes.

The ULTRA X 052 incinerator comes with an electronic with LCD display and 4 switches for the menu and the times and power settings.

The heating time is shown in the display and stage 2 follows automatically when stage 1 is ready. At the end of the process the heater is switched off.

A light and a signal shows when the device is in use.

Use

This infrared rapid incinerator is used particularly in conjunction with ULTRA X laboratory devices to determine the incineration residue in sewage sludge after the determination of dry matter. The UX 052 is useful in every laboratory where a fast incineration of single samples of organic materials is needed.

To determine the dry matter and residue on ignition in sewage sludges, we offer the moisture tester UX 3011 G in combination with the incinerator UX 052.

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Note:

Because noxious and/or malodorous fumes are likely to be emitted during the calcination or ashing process, we recommend that you use the ULTRA X rapid incinerator in a protection cabinet.

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Technical data of the ULTRA X 052

Power supply	230 Volt, 50-60 Hz
Other voltages with separate transformer	
Radiator power:	375 Watt
Temperature range	up to approx. 610 °C
Temperature setting	electronically, 2 steps possible
Time setting	electronically, 2 steps possible up to 99 min. per stage
Sample volume	max. 130 cm ³ (aluminium pan)
Max. bowl size	75 mm diameter x 30 mm height
Dimensions	Height 180 mm Width 190 mm Depth 350 mm
Weight	approx. 4.0 kg

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Accessories available upon request:

Item	Order No.
Aluminium foils 0,03, Ø 130mm	1000017
Foil press to shape the aluminium foils	1000062

Our product range:

Infrared moisture analysers for different usage, sample dryers (without balance), incinerators.